Meditech Europe

MD4 water distiller

user manual

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Unpack the MD4 water distiller and check that all parts are present and undamaged. The package contains:

- Distiller
- Waterjug
- Lid + drain cap
- Filter holder
- Power cable
- Instruction manual
- Carbon filter bag


Assembling:

Our website features an extensive video which explains how to use the MD4

1. Assemble glass water jug

2. Stainless steel water outlet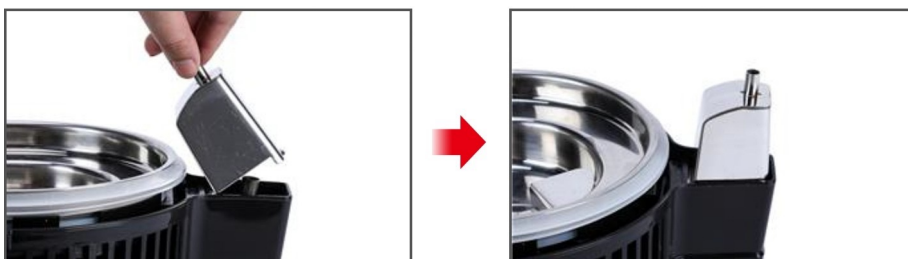

Before use:

Let the MD4 succeed 1 full distillation cycle before using the water.

Cleaning:

1. Rinse and wipe the tank with clean water every day. Calcification and scaling will settle at the bottom of the tank and should be removed after each cycle. Use the following instructions:
2. Pour hot water into the tank until the scale is covered
3. Pour 1-4 teaspoons of vinegar into the tank, depending on the amount of scale.
4. Remove the lid from the MD4 and start a cycle.
5. Let the cycle run for 30 minutes, pour out the water and wash the inside of the tank with clean water.
6. If the scale is not completely removed, repeat the process.
7. Replace the carbon filter every 3 months.

Operation:

1. Fill the tank inside the machine to the highest water level line. Take care not to overfill the tank. (The time needed to distil water can be reduced by using already boiled water).
2. Place the lid on the body of the machine and make sure the silicone seal around the entire edge is firmly in place. Check that the machine is on a stable surface. Then connect the power cable from the top to the MD4 casing.
3. Assemble the waterjug according to diagram I. Connect the power cable to the socket. The fan will turn on automatically and the tank will start heating, starting the distillation process.
4. After about 30 minutes, water will start dripping into the water jug. It takes about 4 hours to distil 4 litres of water. The unit will automatically switch off when the cycle is complete.

Caution:

- Only use the MD4 in a dry, well-ventilated room.
- Keep the outside of the MD4 dry - only fill with water using a jug.
- As soon as the plug is plugged in, the distilling cycle will start. The MD4 will automatically switch itself off when the cycle is complete.
- To start a new cycle, press the reset button This can be found on the side of the MD4.
- Let the MD4 cool down completely before using it again.
- Unplug the MD4 when it is not in use.
- Do not remove the lid when the unit is halfway through its cycle. Wait for the unit to cool down before removing the lid.
- The MD4 must cool down before starting a new distillation cycle. (This prevents damage and prolongs the life of your MD4).

Meditech Europe

Daalder 14

8305 BE Emmeloord

+31 527 292331

info@meditecheurope.nl

www.meditecheurope.nl